


MiniSeq System

Custom Primers Guide

For Research Use Only. Not for use in diagnostic procedures.

Introduction	3
Primer Positions on the Reagent Cartridge	4
Compatible Primer Combinations	5
Prepare and Add Custom Primers	6
Set Up a Run for Custom Primers	7
Revision History	8
Technical Assistance	9

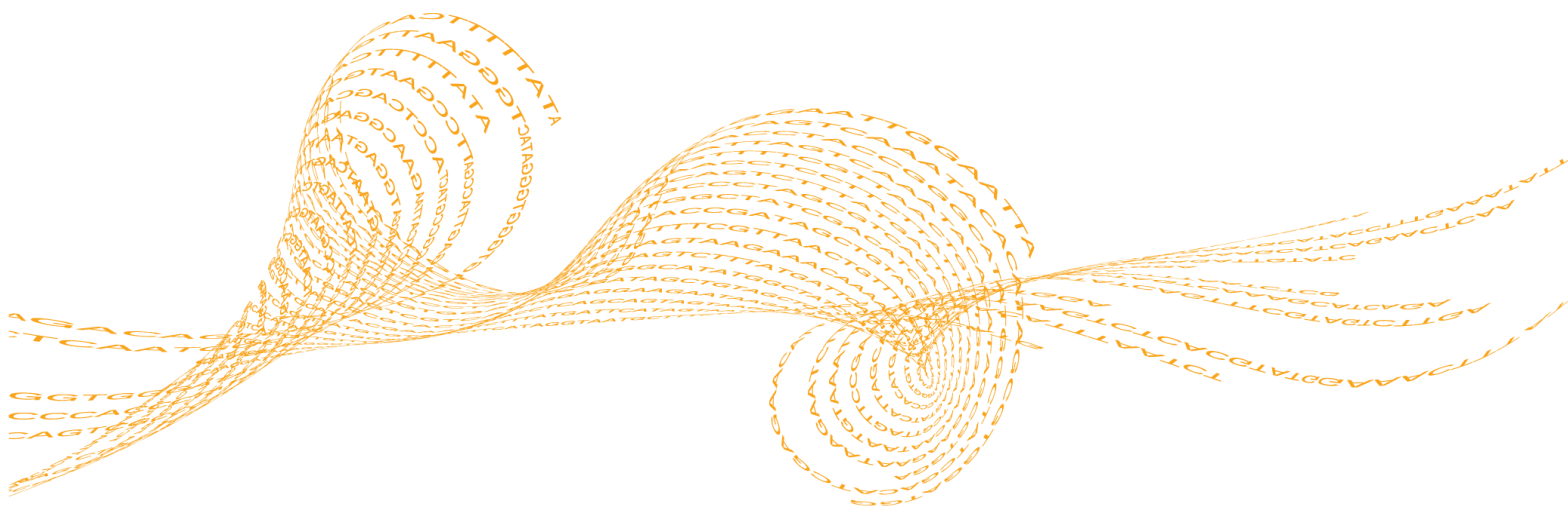

ILLUMINA PROPRIETARY

illumina®

This document and its contents are proprietary to Illumina, Inc. and its affiliates ("Illumina"), and are intended solely for the contractual use of its customer in connection with the use of the product(s) described herein and for no other purpose. This document and its contents shall not be used or distributed for any other purpose and/or otherwise communicated, disclosed, or reproduced in any way whatsoever without the prior written consent of Illumina. Illumina does not convey any license under its patent, trademark, copyright, or common-law rights nor similar rights of any third parties by this document.

The instructions in this document must be strictly and explicitly followed by qualified and properly trained personnel in order to ensure the proper and safe use of the product(s) described herein. All of the contents of this document must be fully read and understood prior to using such product(s).

FAILURE TO COMPLETELY READ AND EXPLICITLY FOLLOW ALL OF THE INSTRUCTIONS CONTAINED HEREIN MAY RESULT IN DAMAGE TO THE PRODUCT(S), INJURY TO PERSONS, INCLUDING TO USERS OR OTHERS, AND DAMAGE TO OTHER PROPERTY.

ILLUMINA DOES NOT ASSUME ANY LIABILITY ARISING OUT OF THE IMPROPER USE OF THE PRODUCT(S) DESCRIBED HEREIN (INCLUDING PARTS THEREOF OR SOFTWARE).

© 2016 Illumina, Inc. All rights reserved.

Illumina, BaseSpace, MiniSeq, the pumpkin orange color, and the streaming bases design are trademarks of Illumina, Inc. and/or its affiliate(s) in the U.S. and/or other countries. All other names, logos, and other trademarks are the property of their respective owners.

Introduction

Using custom sequencing primers for a run on the MiniSeq system requires 2 additional steps during run setup:

- 1 Specify that custom primers are part of the run.
 - ▶ For runs connected to Local Run Manager, use of custom primers are specified on the Create Runs screen.
 - ▶ For runs connected to BaseSpace or BaseSpace Onsite, use of custom primers are specified on the Planned Runs screen of the Prep tab.
 - ▶ For runs using the standalone configuration, use of custom primers are specified on the MiniSeq Control Software Run Setup screen.
- 2 Prepare and add 2 ml of each custom primer to the custom primer positions on the MiniSeq reagent cartridge. See *Prepare Custom Primers* on page 6.

All other steps follow the run setup workflow as described in the *MiniSeq System Guide* (document # 1000000002695).

Custom Primers and PhiX

When custom primers are used for Read 1 or Read 2, the software directs the instrument to pull from reservoirs 14 and 15. Therefore, Illumina primers are not used for the sequencing run.

If Illumina primers are not used for Read 1 or Read 2, the optional Illumina PhiX control is *not* sequenced. To use the PhiX control with custom primers, contact Illumina Technical Support for guidance.



NOTE

Sequencing data from the PhiX control is not generated for index reads, regardless of which indexing primer is used.

Primer Positions on the Reagent Cartridge

You can use a combination of Illumina primers and custom primers in the same run. Depending on the combination specified, the software pulls the primer from the appropriate reservoir.

Example: If a custom primer is used for Read 2, but not for Read 1, the software pulls the Read 1 primer from Illumina Read 1 primer position (#24) and pulls the Read 2 primer from the custom Read 2 primer position (#14).

Custom Primer	Custom Primer Position
Custom Read 1 primer	15
Custom Read 2 primer	14
Custom Index 1 primer	13
Custom Index 2 primer	

Illumina Primer	Illumina Primer Position
Illumina Read 1 primer	24
Illumina Read 2 primer	25
Illumina Index 1 primer and Index 2 primer mix	28

Primer positions used for the run are recorded in the file Primers.xml, which is written to the Reagents folder.

Compatible Primer Combinations

Due to the primer design and dual-indexing workflow on the MiniSeq, combinations of custom primers and Illumina primers are limited.

- ▶ If you use a custom Read 1 primer, you have to use a custom Index 2 primer.
- ▶ If you use a custom Read 2 primer, you have to use a custom Index 1 primer.

For an overview of the indexing workflow, see *Indexed Sequencing Overview Guide* (document # 15057455).

Combination	Read 1 Primer	Index 1 Primer	Index 2 Primer	Read 2 Primer
All Illumina	Illumina Read 1	Illumina Index 1	Illumina Index 2	Illumina Read 2
Custom Read 1	Custom Read 1	Illumina Index 1	Custom Index 2	Illumina Read 2
Custom Read 2	Illumina Read 1	Custom Index 1	Illumina Index 2	Custom Read 2
All custom	Custom Read 1	Custom Index 1	Custom Index 2	Custom Read 2

Prepare and Add Custom Primers

Make sure that the MiniSeq reagent cartridge is thawed and inspected before proceeding.

Prepare Custom Primers

- 1 If frozen, thaw each custom primer to be used.
- 2 Use HT1 to dilute each custom primer to yield 2 ml at 0.3 μ M final concentration.

Add Custom Primers to Reagent Cartridge

- 1 Use a low-lint lab tissue to wipe clean the foil seal covering each position associated with the custom primer.

Position #	Custom Primer
15	Custom Read 1 primer
14	Custom Read 2 primer
13	Custom Index 1 primer Custom Index 2 primer

- 2 With a clean pipette tip, pierce the foil seal covering each position associated with the custom primer.
- 3 Add the following volumes of custom primer in the appropriate position on the reagent cartridge. Avoid touching the foil seal as you dispense the primer.

Position #	Volume	Custom Primer
15	1 ml	Custom Read 1 primer
14	1 ml	Custom Read 2 primer
13	1 ml	Custom Index 1 primer
13	0.5 ml of each primer	Custom Index 1 primer <i>and</i> Custom Index 2 primer
13	1 ml	Custom Index 2 primer

Set Up a Run for Custom Primers

The use of custom primers is specified during run setup.

- ▶ If the instrument is configured for Local Run Manager, specify custom primers on the Create Run page.
- ▶ If the instrument is configured for BaseSpace or BaseSpace Onsite, specify custom primers on the BaseSpace Prep tab. You can later edit these settings on the MiniSeq Control Software Run Setup screen.
- ▶ If the instrument is configured for standalone mode, specify custom primers on the MiniSeq Control Software Run Setup screen.

Specify Custom Primers Using BaseSpace Prep Tab

- 1 On the Plan Run screen of the BaseSpace Prep tab, select the checkbox for each custom primer to be used with the run.
 - ▶ **R1**—Custom primer for Read 1.
 - ▶ **R2**—Custom primer for Read 2.
 - ▶ **Index**—Custom primer for Index 1 Read.To specify custom primers for Index 1 or Index 2 separately, edit the custom primers settings on the MiniSeq Control Software Run Setup screen.

Specify Custom Primers Using Local Run Manager

- 1 On the Create Run page of Local Run Manager, select the checkbox for each custom primer to be used with the run.
 - ▶ **Read 1**—Custom primer for Read 1.
 - ▶ **Index 1**—Custom primer for Index 1 Read.
 - ▶ **Index 2**—Custom primer for Index 2 Read.
 - ▶ **Read 2**—Custom primer for Read 2.

Specify Custom Primers in Standalone Configuration

- 1 On the MiniSeq Control Software Run Setup screen, select the checkbox for each custom primer to be used with the run.

During the run, the software directs the sippers to pull primers from the appropriate reservoir.
- 2 When you have finished specifying run parameters, select **Next**.

Revision History

Document	Date	Description of Change
Document # 1000000002700 v01	September 2016	Updated custom primer positions to 15 for Read 1, 14 for Read 2, and 13 for Index 1 and Index 2.
Document # 1000000002700 v00	February 2016	Initial release.

Technical Assistance

For technical assistance, contact Illumina Technical Support.

Table 1 Illumina General Contact Information

Website	www.illumina.com
Email	techsupport@illumina.com

Table 2 Illumina Customer Support Telephone Numbers

Region	Contact Number	Region	Contact Number
North America	1.800.809.4566	Japan	0800.111.5011
Australia	1.800.775.688	Netherlands	0800.0223859
Austria	0800.296575	New Zealand	0800.451.650
Belgium	0800.81102	Norway	800.16836
China	400.635.9898	Singapore	1.800.579.2745
Denmark	80882346	Spain	900.812168
Finland	0800.918363	Sweden	020790181
France	0800.911850	Switzerland	0800.563118
Germany	0800.180.8994	Taiwan	00806651752
Hong Kong	800960230	United Kingdom	0800.917.0041
Ireland	1.800.812949	Other countries	+44.1799.534000
Italy	800.874909		

Safety data sheets (SDSs)—Available on the Illumina website at support.illumina.com/sds.html.

Product documentation—Available for download in PDF from the Illumina website. Go to support.illumina.com, select a product, then select **Documentation & Literature**.



Illumina

5200 Illumina Way

San Diego, California 92122 U.S.A.

+1.800.809.ILMN (4566)

+1.858.202.4566 (outside North America)

techsupport@illumina.com

www.illumina.com